


Automatic Leveling Bracket Reinstallation

1. Turn the Onyx table upside down, exposing the underside of the table.
2. Remove all parts of the old blue bracket that remains on the valve by unbolting the blue bracket from the existing valve.
3. Remove the pin from the lever arm.
4. Insert the lever arm into the new bracket as shown in the diagram below.
5. Place the pin into the bracket and lever arm holes.
6. Attach the bracket assembly to the valve.
7. Adjust level of the system per the instructions found in the Onyx Series Automatic Leveling Manual.

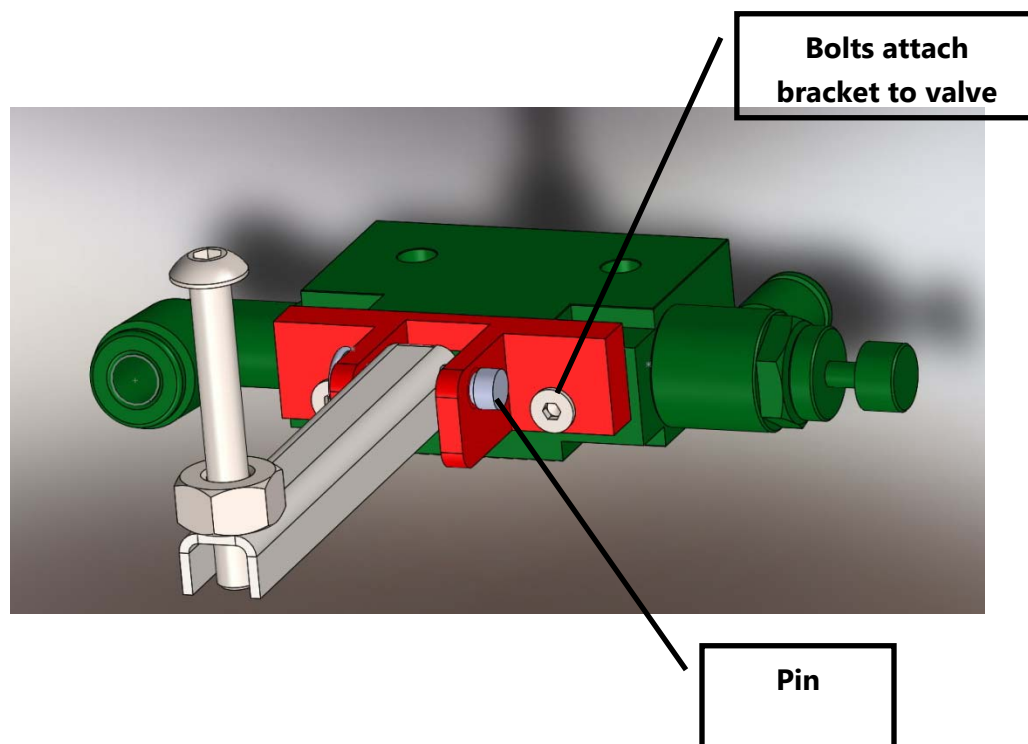

Troubleshooting

For troubleshooting information: <http://herzan.com/support/troubleshooting/air-based-isolation.html> or follow the QR code on your mobile device provided below:



Any Questions?

Contact us today if you have any questions or would like assistance with your installation. Herzan is available Monday – Friday, from 8:00 A.M. to 6:00 P.M. (Pacific Standard Time).

T: (949) 363-2905

E: support@herzan.com

